

Industrial Legionella TestTM

...delivering microbiology results in minutes instead of days

- Fast results
- High sensitivity
- In-line sampling
- Ideal for industrial samples

The Industrial Legionella TestTM, is a breakthrough technology from hydrosense[®] for the rapid detection of *Legionella* bacteria in industrial water – delivering test results in minutes instead of days.

The test kit contains all the items required to perform 5 tests – including filters, recovery buffer, collection vessel and sample point.

- These tests are optimised for use with **industrial** water systems, **cooling towers** and whirlpool **spas**.
- This filtration step needs **no pumping**, just connect the sample point to the water system and use the system pressure to drive the filtration.
- Single use **hollow fibre filtration modules**.
- Pre-filled syringe containing recovery buffer, no need to measure or dilute solutions.
- The test is **simple to use** and **fast**, with no interpretation required.
- Results are **easy to read** and obtained within 30 to 35 minutes.
- The test allows **immediate understanding** and response to the water conditions.
- Use can **reduce risk** of outbreaks and litigation.
- Regular use will **increase confidence** in your *Legionella* control.



Industrial Legionella TestTM
Product code 100182

The hydrosense[®] Legionella Field TestTM uses an immunochromatographic assay, to detect the presence of cell surface antigens from *Legionella* bacteria in a sample. The presence of antigen causes the “Test Line” to turn red in colour. A “Control Line” is included which should always turn red if the test has been performed correctly.

Product details

Bacteria detected *Legionella pneumophila* serogroup 1

Detection method Lateral flow immunochromatographic assay

Lower limit of detection Up to 100 CFU/L (0.1 CFU/mL)

Specificity

Tested against a wide range of waterborne bacteria, including *Acinetobacter calcoaceticus*, *Aeromonas hydrophila subsp. hydrophila*, *Bacillus subtilis*, *Burkholderia cepacia*, *Citrobacter freundii*, *Citrobacter koseri*, *Escherichia coli*, *Enterobacter cloacae*, *Klebsiella oxytoca*, *Pseudomonas aeruginosa*, *Pseudomonas fluorescens*, *Pseudomonas putida*, *Pseudomonas stutzeri*, *Ralstonia pickettii*, *Raoultella terrigena*, *Streptococcus pyrogenes*, *Yersinia ruckeri*. No cross reactions observed at concentrations $\leq 10^9$ CFU/L. *Staphylococcus aureus* has been observed to cross react with the test at concentrations $> 10^8$ CFU/L. *Legionella pneumophila* serogroups 4 & 7 cross react with the test at concentrations $> 10^9$ CFU/sample. The test does not react with any other *L. pneumophila* serogroups or *L. species* tested.

Test volume 100 μ L

Suggested sample volume 250 mL

Sample parameters Sample should filter in less than 10 minutes.

Biocides


The product has been tested with a range of common biocides and biodispersants used in *Legionella* control, at their normal maximum operating concentration. Active ingredients tested include: glutaraldehyde, didecyl-dimethyl-ammonium chloride, isothiazoles, dibromonitropropionamide, ionic and non-ionic surfactants. The test gives a false positive if used with polymeric biguanide or THPS.

Operator time 5 - 10 minutes

Test time 25 minutes

Sample point Supplied with both 1/2" and 3/4" female threaded fittings, additional sample points are available (part no. 100171)

Shelf life 18 months from manufacture when stored at room temperature ($< 30^{\circ}\text{C}$) in original packaging.

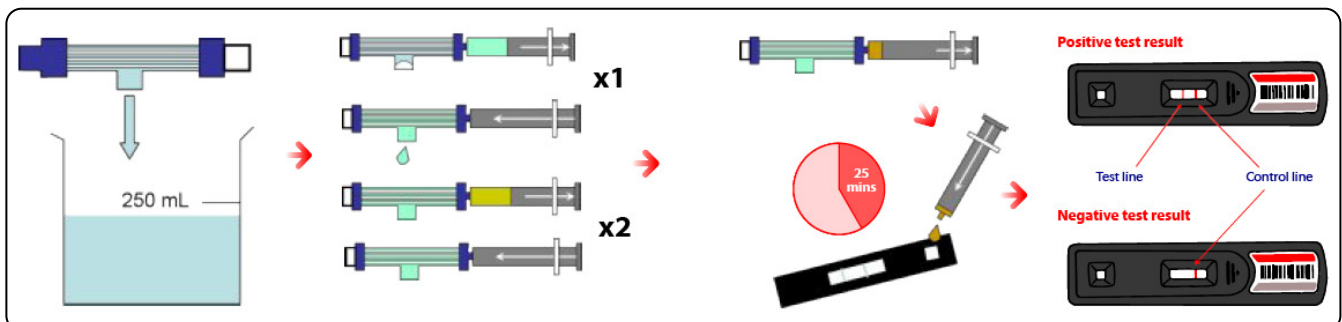

Albagaia Limited Pinnacle House Linlithgow EH49 7SF United Kingdom
Phone: +44 (0)1506 841804 Fax +44 (0) 871 2668162

www.hydrosense.biz

TECH0019v0.1

hydrosense®